

Oscilloscope Set-up For Wohlwend Compact 02 High Pressure Freezer

Software Installation and Basic set-up for Windows 7

First time only:

Connect PC to USB port via enclosed USB cable


Start Wohlwend Compact 02 HPF

Start-up PC

Windows will discover new devices, one for channel A and one for channel B.

Then install EasyScope software from supplied USB stick

Start-up EasyScope from desktop icon

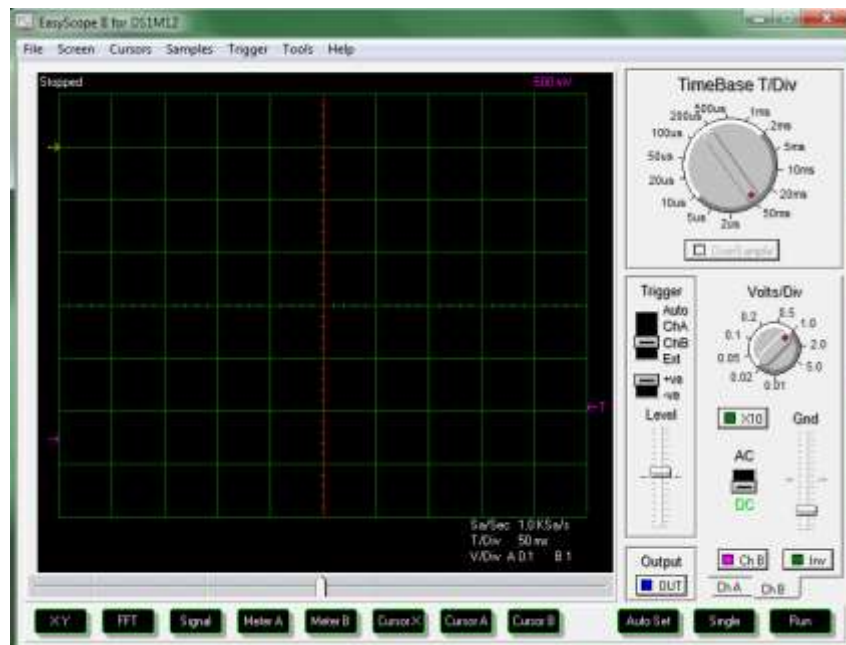

Oscilloscope Settings:

TimeBase T/Div:	50 ms	
Channel A (ChA):	0.1 Volts/Div	measures temperature default line yellow
Channel B (ChB):	0.1 Volts/Div	measures pressure default line color purple

To set the trigger, slide Trigger switch to ChB, and set trigger point to 600 mV

Operation:

Push single on oscilloscope, the screen will go blank
Push Jet Auto on System

With the slide below the screen, the curve can be shifted from left to right

Save the screen image:

Under the screen menu, save screen image, the curve can be saved either into the default directory or a directory of your choice.

Default directory:

C://ProgramFiles/USBInstruments/Easyscope/Images

Link to manufacture's website for documentation and software / driver downloads:

<http://www.easysync-ltd.com/product/520/ds1m12.html>

TECHNOTRADE International, Inc. | 7 Perimeter Road |
Manchester, NH | 03103 (603) 622-5011
sales@technotradinc.com