

PROD. NO. 607
MAGNIFICATION CALIBRATION
DIFFRACTION GRATING REPLICA

CAT # 603

Latex Spheres on Diffraction Grating Replica

This specimen is a replica of a 2,160 lines/mm waffle-pattern diffraction grating on which 0.261µm diameter latex particles have been applied. When imaging the specimen it should be kept in mind that the line spacing is 0.463µm and the pattern will not be visible until the imaging system is set to resolve that level of detail - around x2,500. At this magnification the lines of the pattern will be 1.1575 mm apart and the latex particles will have a diameter of .625 mm.


1. To calculate the electron microscope magnification using the pattern of the **diffraction grating replica**:

Take the measurement, in millimeters, between the limiting lines of as many squares of the replica pattern as possible. Apply the following formula:

$$\text{Magnification} = A \times 2,160/B$$

A is distance in mm between limiting lines.
B is number of spaces between limiting lines.

[Alternatively, use the PelcoTM Magnification Calibration Calculator, Cat. No. 252]

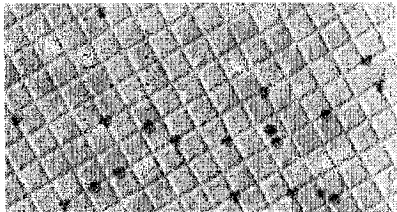

2. To calculate the magnification by measuring the magnified image of the 0.262µm dia. latex spheres, apply the following formula:

$$\text{Magnification} = \text{diameter (in mm)} \times 3,831.4$$

Note: Due to variations in size, the latex particles are not an accurate way of determining instrument magnification, although the calculated figure will be tolerably close to that obtained using the diffraction grating pattern. They serve rather, as a useful point of reference for visualizing the appearance of objects at different magnifications and determining the lowest magnification at which structures you are looking for, in another specimen, might be visible.

Care of Grating Replica Specimen

When not in use the replica should be kept in the vial. The replica surface may be damaged if touched. Never try to clean it. Care must be taken to avoid bending the grid as distortion may cause the replica film to fracture. **NOTE: ALWAYS VIEW THIS SPECIMEN AT THE LOWEST COMFORTABLE ILLUMINATION LEVEL TO AVOID DEGRADATION OF THE LATEX PARTICLES.**




Ted Pella Inc.

Tools for Science

P.O. Box 492477, Redding California 96049-2477

Phone: (530) 243-2200 Fax: (530) 243-3761

This specimen is a replica of a 2,160 lines/mm waffle-pattern diffraction grating. When imaging the specimen it should be kept in mind that the line spacing is 0.463µm; also the pattern will not be visible until the imaging system is set to resolve that level of detail, which is around x2,500. At this magnification the lines of the pattern will be just over 1mm apart.

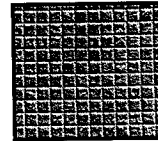

To calculate the electron microscope magnification using the pattern of the diffraction grating replica:

Take the measurement, in millimeters, between the limiting lines of as many squares of the replica pattern as possible. Apply the following formula:

$$\text{Magnification} = A \times 2160/B$$

A is distance in mm between limiting lines.
B is number of spaces between limiting lines.
(Alternatively, use the Pelco[®] Magnification Calibration Calculator, Prod. No. 252.)

Care of Grating Replica Specimen

When not in use, the replica should be kept in the vial. The replica surface may be damaged if touched. Never try to clean it. Care must be taken to avoid bending the grid as distortion may cause the replica film to fracture. When viewing in the TEM begin at low magnification with a low illumination level. Increase the illumination a little beyond comfortable viewing level then reduce it. This helps to stabilize the specimen. Before moving the specimen to view another grid square, reduce the illumination and magnification to starting levels again.

Note:

Artifacts such as furrows or other distortions are on the gold surface of the original master grating and do not affect the accuracy of the line spacing

607 TN V3 06032005

TED PELLA, INC.
Microscopy Products for Science and Industry

PO Box 492477, Redding, CA 96049-2477
Fax (530) 243-3761 Phone: (530) 243-2200
(800) 237-3526 (USA and Canada)

Email: sales@tedpella.com Website: www.tedpella.com